

protocol_22

Table of contents

1	Protocol Parameters	1
2	Conditions	2
3	Pattern Previews	2
3.0.1	Pattern 31	2
3.0.2	Pattern 32	2
3.0.3	Pattern 33	2
3.0.4	Pattern 34	3
3.0.5	Pattern 35	3
3.0.6	Pattern 36	3
3.0.7	Pattern 37	4
3.0.8	Pattern 38	4

Protocol_22.m - follow up protocol 2 after protocol 19 double effector test screen. Based on protocol 19 Includes: - greyscale interval frame. - 4 pixel OFF bar fixation - 180 deg apart - Reverse Phi - 4 bit - 1 pix step - 8 pixel width bar - 24 px gap. - Reverse Phi - 4 bit - 4 pix step - Reverse Phi - 4 bit - 8 pix step

1 Protocol Parameters

Parameter	Value
Conditions	8
Patterns Used	8
t_acclim	20s
t_pause	0.015

2 Conditions

#	Pattern	Speed	Duration	Description
1	Pattern 37	29	-s	1 = bar fixation - OFF bar
2	Pattern 38	29	-s	1 = bar fixation - ON bar
3	Pattern 31	29	-s	reverse phi - 1px step
4	Pattern 32	29	-s	4 px step
5	Pattern 33	29	-s	8 px step
6	Pattern 34	29	-s	FoE - 30deg
7	Pattern 35	29	-s	15 deg
8	Pattern 36	29	-s	60 deg

3 Pattern Previews

3.0.1 [Pattern 31](#)



3.0.2 [Pattern 32](#)




3.0.3 [Pattern 33](#)




3.0.4 Pattern 34



3.0.5 Pattern 35

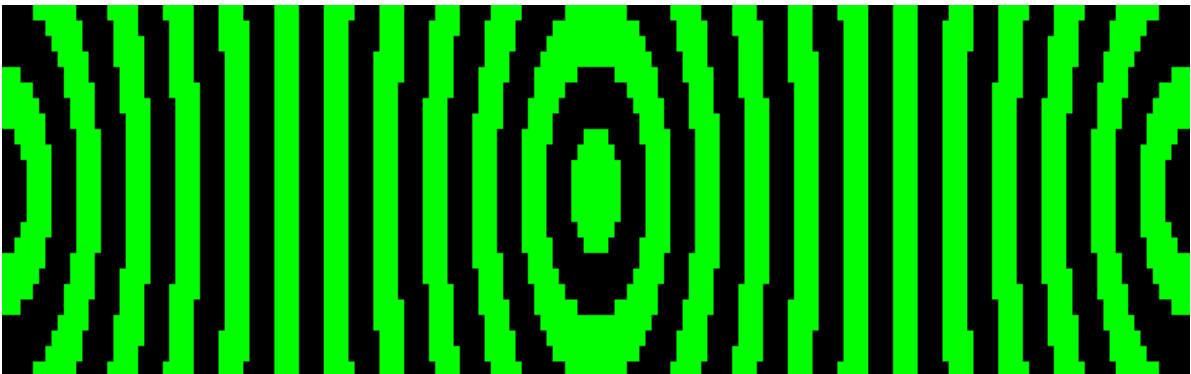

3.0.6 Pattern 36

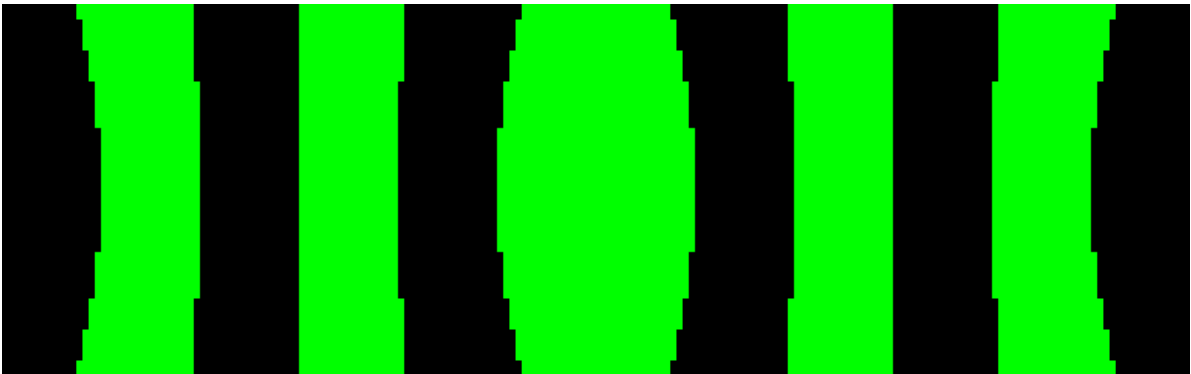

3.0.7 **Pattern 37**



3.0.8 **Pattern 38**

