

protocol_26

Table of contents

1	Protocol Parameters	1
2	Conditions	2
3	Pattern Previews	2
3.0.1	Pattern 47	2
3.0.2	Pattern 51	2
3.0.3	Pattern 52	2
3.0.4	Pattern 55	2
3.0.5	Pattern 56	2

Protocol_26.m - test different curtain stimuli. 32 pixels instead of 16. cond 1 = binary on cond 2 = binary off cond 3 = greyscale ON cond 4 = greyscale OFF 20s interval 30s trial duration - 2x trials

1 Protocol Parameters

Parameter	Value
Conditions	4
Patterns Used	6
t_acclim_start	30s
t_acclim_end	30s
t_interval	20s
t_flash	5s
t_pause	0.01

2 Conditions

#	Pattern	Speed	Duration	Description
1	Pattern 51	127	15s	ON - 32px - binary - 1-0
2	Pattern 52	127	15s	OFF - 32 px - binary - 1-0
3	Pattern 56	127	15s	ON - 32px - greyscale - 4-0
4	Pattern 55	127	15s	OFF - 32px - greyscale - 4-0

3 Pattern Previews

3.0.1 [Pattern 47](#)



3.0.2 [Pattern 51](#)



3.0.3 [Pattern 52](#)



3.0.4 [Pattern 55](#)



3.0.5 [Pattern 56](#)

